

## ○ Virus

## ● Marburg virus

The NP gene of Marburg virus, the causative agent of Marburg virus disease, was detected by ultra-rapid real-time RT-PCR.

## ○ Instruments and reagents

- GeneSoC® mini R
- GeneSoC® mini PCR chips
- GeneSoC® PCR RT-Master Mix

## ○ Primers and Probes


J Clin Lab Anal. 2021;35(6):e23786.

## ○ Template

Synthetic RNA (Marburg virus NP gene)

## Related Products

- Ultra-rapid real-time PCR system GeneSoC® mini R
- GeneSoC® mini PCR chips
- GeneSoC® RT-PCR Master Mix

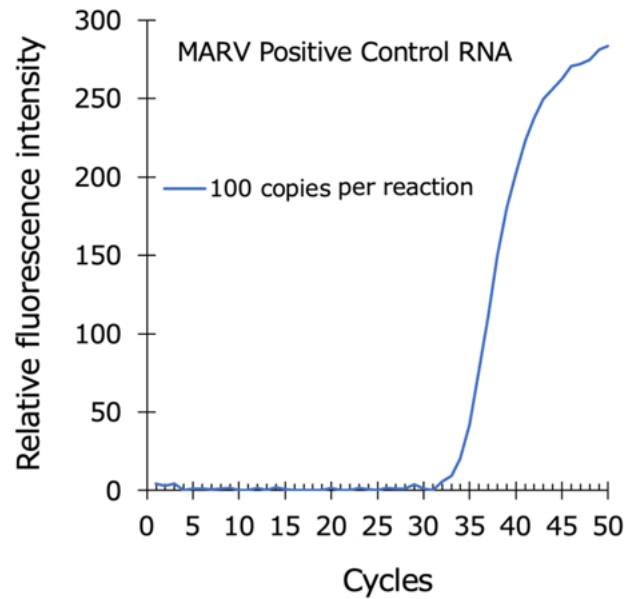

## ○ PCR Protocol

	Temp.(°C)	Time(sec)	Cycle
RT	42	60	1
HS	98	5	
DN	98	3	50
AE	60	8	

(Time required: approx. 14.5 minutes)

- HS : Hot Start
- DN : Denature
- AE : Annealing / Extension

※ For more detailed experimental conditions, please contact us.



Ultra-rapid real-time PCR system

**GeneSoC® mini R**

<https://genesoc.jp/en/application-note/>



Palm-sized PCR device

